

BIODENE High performance Hydrophilic PVDF

- Easy integrity testable in situ
- Repeatedly steamable in situ and in autoclave
- Thermowelded construction
- EC-listed materials for Food contact
- FDA-listed materials per 21 CFR
- Bio-Safety per USP—Plastics



The new BIODENE filter element is a perfect combination between the PVDF membrane with enhanced hydrophilic characteristic and the polypropylene stratification with its high mechanical resistance. The BIODENE filter element has been designed to meet the high standard quality and safety requirements for the most critical application of oenology and beverages applications.

This filter element assures the microbiology stability and remove all the contaminants not welcome during the bottling activity, keeping the original color, flavor and scent of the filtered product.

The controlled porosity together with the high resistance to heat sterilization, the chemical compatibility and the resistance to mechanical strengths are the highlights of this new filter elements series.

Its enhanced service life and the easy regenerations activity even using soda, contributes to a great cost reduction specially when treating fluids with high colloidal burden.

Typical applications are all the soft drinks, water, wine, cider and raw beer.

The manufacturing is performed in a controlled environment; all the filter elements are 100% integrity tested and verified in production.

The filter element is available in 0.2, 0.45 and 0.65 micron configuration.

MATERIALS OF CONSTRUCTION

Filter media	Hydrophilic PVDF	
Upstream supports	polypropylene	
Downstream supports	polypropylene	
Internal Core	polypropylene	
External Cage	polypropylene	
End caps	polypropylene	

FOOD-SAFETY

BIODENE filter element materials meet (EU) regulation 10/2011 and its amendments, regulations (EC) 1935/2004 and 1895/2005.

BIO-SAFETY

Filter media and components pass USP CLASS VI Biological Reactivity and Chemical-Physical tests for USP plastics.

QUALITY STANDARDS

Produced under a certified Quality System to guarantee traceability of manufacturing records and integrity testing results.

RECOMMENDED OPERATING CONDITIONS

- max. continuous temperature	85 °C
- max. cumulative time of steam sterilization	80 hours at 125 °C with cycles of 60 minutes />100 hours at 121 °C
- sanitization with hot water	90 °C max
- sanitization with chemicals	Can be sanitized by standard chemical agents
- max. differential pressure	5,0 bar at 25 °C—2,5 bar 80 °C
- recommended change out differential pressure	2,0 bar at 25 °C
- regeneration	NaOH solution up to 2% at 80 °C

CODE	CODE ABSOLUTE FILTRATION RATING IN LIQUIDS	MAX. DECAY VALUE *	ACCEPTABLE LIMIT FOR DIFFUSION FLOW TEST WITH WATER FOR 10" CARTRIDGE
		8 of 30" CARTRIDGES	(ml/min)
BSY	0,2 μm	≤ 0,12 bar	≤ 25 @ 1,6 bar
BST	0,45 μm	≤ 0,12 bar	≤ 25 @ 1,2 bar
BSK	0,65 μm	≤ 0,11 bar	≤ 22 @ 0,9 bar
BSF	1,0 μm	≤ 0,10 bar	≤ 20 @ 0,5 bar

The values are related to 5 minutes and are indicative as they depend on the housing volume upstream the filter element.

MAX FLOW RATE FOR 30" CARTRIDGE


CODE	L/H
BSY	1000
BST	1500
BSK	2000
BSF	2500

BACTERIAL RETENTION


CODE	FILTRATION RATING	*BACTERIAL RETENTION > 10 ⁷ per cm ²
BSY	0,2 μm	Pseudomonas aeruginosa / Escherichia coli / Enterobacteriacee
BST	0,45 μm	Oenococcus oeni/ Saccharomyces cerevisiae / Brettanomyces bruxellensis / Lactobacilius brevis / Oocystes cryptosporidium / Giardia / Pediococcus damnosus
BSK	0,65 μm	Saccharomyces cerevisiae/ Brettanomyces bruxellensis
BSF	1,0 µm	Saccharomyces cerevisiae/ Brettanomyces bruxellensis (10 ⁶ per cm ²)

^{*} Secondo ASTM F838

WATER FLOW RATE CURVES FOR 10" ELEMENT

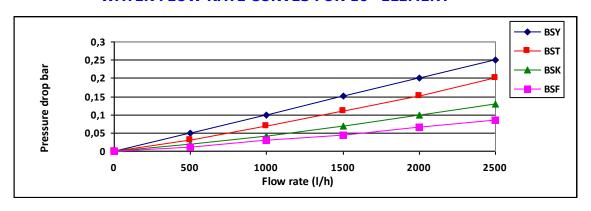

BIODENE ORDERING INFORMATION

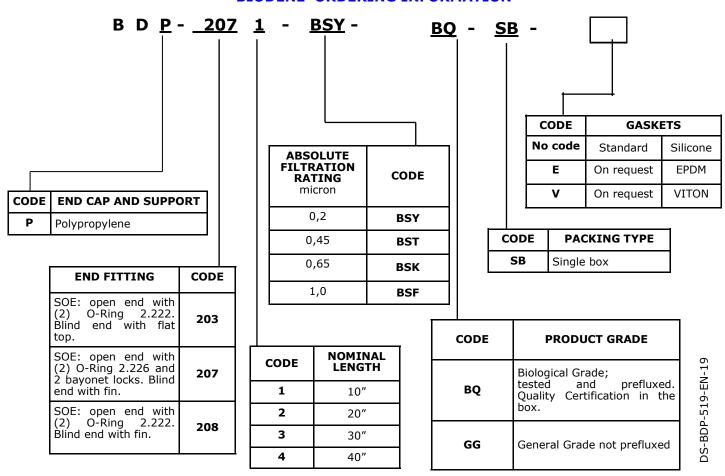

Data contained in this bulletin are informative and subject to change without notice. User is responsible for determining whether the product is fit for particular purpose and suitable for User's method of application.



Bea Technologies Spa Via Newton, 4 - 20016 Pero (Milano) ITALY Tel +39 02 339271 FAX +39 02 3390713 e-mail: info@bea-italy.com web: www.bea-italy.com